

## LOTF Laser Shutdown Procedure

### S-AB-P-039 Rev A

**Precursor:**

This is a Class 3B laser operating at 1053nm. Appropriate eye protection should be worn.

**Shutdown Procedure:**

- 1 Turn power knob fully counter-clockwise until it stops and power reads 0 mW.

---

  - 2 Turn key on rear of unit to its 'off' position.

---

  - 3 Inform the Shot Director that the LOTF laser has been turned off.
- 

(End of procedure)

**S-AB-P-039 LOTF Laser Startup Procedure**

**Document Release:**

This document is a component of Vol. IX OMEGA EP System Startup and Shut Down, Chapter 4, Beamlines Operating Procedures, S-AB-P-013.

Approval for release of this document into the PDM system was granted by:

S. Loucks; Director, LLE Engineering Division  
S. Morse; OMEGA EP Project Manager  
J. Edwards; OMEGA EP System Engineer  
R. Jungquist; OMEGA EP Beamlines Subsystem Engineer  
M. Moore; Beamlines Engineer, LOTF  
A. Rigatti; Manager, Optical Manufacturing